

Ultraclean SA Technical Data Sheet

Description:

Ultraclean SA is a mild detergent for general use in ultrasonic baths.

- Aqueous based product
- General purpose cleaning
- Reduces surface tension, enhances ultrasonic cleaning performance
- Contains de-foaming agent
- Suitable for use on aluminium

Applications

Used in a wide range of applications including instrument decontamination in healthcare and cleaning different components in manufacturing and processing industries. Ultraclean SA can be used to remove contaminants from plastics, glass, metals and other materials.

Directions for Use

Add 20-50ml of Ultraclean SA per 1 Litre of water in the ultrasonic cleaner. Do not mix with other solutions.

The required dosage will depend on your cleaning application and the level of contamination. Items should be rinsed with water and dried immediately after processing.

Technical Data

These figures are typical and should not be interpreted as specifications, refer to MSDS for details.

Characteristic	Value/Description
Appearance	Green liquid
Odour	Mild, characteristic
pH (@ 1%)	9.8 (typical)
Density at 20°C	1.112



Available in:

- 6x1L box (Part no. F5002501)
- 25 Litre Drum (Part no. F5002401)

Disposal

When diluted to concentration of 5% or less, Ultraclean SA is not hazardous to aquatic organisms and may be safely disposed of down the drain. Refer to regional or national regulations.

Storage

12 months when stored in a cool, dark environment.